

Capillary Adapter NE-382-CAOR Assembly and Use

Insert the Bushing into the base of the syringe as shown.

The nut holds the O-ring as shown. Make sure that the O-ring is seated properly.

Thread the screw onto the syringe, taking care not to cross-thread it, until it is firmly in place. Do not over tighten.

Taking note of the alignment marks on the screw and nut, thread the nut onto the screw until you feel the slightest resistance. The first line on the nut should align with the line on the screw. This is the open position where the glass needle can be easily inserted in place. The needle will stop at the bushing.

A small turn clockwise would secure the needle. The second mark on the nut should align with the line on the screw. A small turn counter-clockwise would release the needle.

The alignment marks are not permanent. They are there initially to help you get a feel for work with the system. There is no need to fully remove the nut except for replacement or cleaning.

